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**Instructions for use
Eye wash bottles
BARIKOS KS and
Mini-BARIKOS KS**

ENGLISH

The filled eye wash bottles BARIKOS complying with EN 15154-4 are intended as first aid method to rinse the eyes with water in order to remove penetrated substances before starting medical treatment.

If required

- **remove the top**
- **bend the head forward over the bottle or put the bottle in front of the injured eye**
- **open the eye with thumb and index finger as far as possible**
- **press the spray-head gently against the eye socket while keeping the eye open and rinse plentifully by squeezing the bottle repeatedly.**

When closing the eyelids reflexly the injured person mostly is not able to rinse by himself sufficiently. In this case the nearest helper has to give first aid.

The eye wash bottle and the intactness of the seal has to be inspected regularly.

After using the bottle the remaining water is to be rejected. The same applies in case the bottle is being opened (broken seal) but not used.

Shelf life of the rinsing liquid (water): 2 years if the bottle remains closed and the seal is unbroken.

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